
Product Manual


OxiSelect™ Protein Carbonyl Fluorometric Assay

Catalog Number

STA-307

100 assays

FOR RESEARCH USE ONLY
Not for use in diagnostic procedures

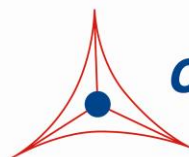

CELL BIOLABS, INC.
Creating Solutions for Life Science Research

Introduction

Protein oxidation is defined as the covalent modification of a protein induced either directly by reactive oxygen species (ROS) or indirectly by reaction with secondary by-products of oxidative stress. Oxidative modification of proteins can be induced in vitro by a wide array of pro-oxidant agents and occurs in vivo during aging and in certain disease conditions.

There are numerous types of protein oxidative modification. The most common products of protein oxidation in biological samples are the protein carbonyl derivatives of Pro, Arg, Lys, and Thr. These derivatives are chemically stable and serve as markers of oxidative stress for most types of ROS. Traditionally, carbonyls have been measured spectrophotometrically through derivatization of the carbonyl group with dinitrophenylhydrazine, although this method has limited sensitivity. In response to this, a new fluorescent compound has been developed this will bind specifically to carbonyls of oxidized proteins at 1:1 molar ratio. Upon binding with the protein carbonyl, the fluorescence produced is a direct measure of protein carbonyls and is more sensitive than the traditional spectrophotometric assay using DNPH.

Cell Biolabs' OxiSelect™ Protein Carbonyl Fluorometric Assay provides a convenient system to measure the protein carbonyl content in plasma, serum, cell lysates, tissues and purified proteins. Each kit provides sufficient reagents to perform up to 100 assays including unknown protein samples, assay blanks and standards.

Assay Principle

Protein carbonyls in protein samples (1-10 mg/mL) are first derivatized with the Protein Carbonyl Fluorophore. The Fluorophore binds to the protein carbonyl group in a 1:1 ratio. Proteins are then TCA precipitated and free Fluorophore is removed by washing the protein pellet with acetone. After dissolving the protein pellet in GuHCl, the absorbance of protein-fluorophore product is measured fluorometrically with a 485/538 nm filter set, and the protein carbonyl is subsequently calculated.

Related Products

1. STA-305: OxiSelect™ Nitrotyrosine ELISA Kit
2. STA-308: OxiSelect™ Protein Carbonyl Immunoblot Kit
3. STA-310: OxiSelect™ Protein Carbonyl ELISA Kit
4. STA-816: OxiSelect™ N-epsilon-(Carboxymethyl) Lysine (CML) Competitive ELISA Kit
5. STA-817: OxiSelect™ Advanced Glycation End Products (AGE) Competitive ELISA Kit

Kit Components

1. Protein Carbonyl Fluorophore (100X) (Part No. 230701): One 50 µL amber vial
2. Sample Diluent (10X) (Part No. 230702): One 20 mL bottle
3. Protein Solubilization Solution (Part No. 231504): One 15 mL bottle
4. Assay Diluent (4X) (Part No. 234505): One 50 mL bottle
5. Fluorophore Standard (Part No. 230703): One 20 µL amber vial of a 20 mM solution in DMSO

Materials Not Supplied

1. 1 to 10 mg/mL protein samples such as plasma, serum, cell lysate, purified protein
2. Protein assay such as Pierce BCA or Bradford assay
3. 20% (w/v) Trichloroacetic Acid Solution (TCA)
4. Acetone
5. Streptomycin sulfate
6. 5 μ L to 1000 μ L adjustable single channel precision micropipettes with disposable tips
7. 50 μ L to 300 μ L adjustable multichannel micropipette with disposable tips
8. Bottles, flasks, and conical or microcentrifuge tubes necessary for reagent preparation
9. Centrifugal filter or concentrator
10. Reagents and materials necessary for sample extraction and purification
11. 96-well black or fluorescence microtiter plate
12. Fluorescent microplate reader capable of reading 480 nm (excitation) and 530 nm (emission)

Storage

Upon receipt, aliquot and store the Protein Carbonyl Fluorophore (100X) and Fluorophore Standard at -20°C to avoid multiple freeze/thaw cycles. Store all other components at 4°C .

Preparation of Reagents

- 1X Sample Diluent: Dilute the provided Sample Diluent (10X) to 1X with deionized water. Stir to homogeneity. Store the 1X Sample Diluent at 4°C .
- 1X Assay Diluent: Dilute the provided Assay Diluent (4X) solution 1:4 with deionized water. Mix to homogeneity. Store the 1X Assay Diluent at 4°C .
- 1X Protein Carbonyl Fluorophore: Prepare only enough for immediate applications. Prepare 1X Protein Carbonyl Fluorophore by diluting the provided Protein Carbonyl Fluorophore (100X) stock solution 1:100 with 1X Sample Diluent. Do not store diluted solution.

Preparation of Samples

All samples should be assayed immediately or stored at -80°C . The assay can be used on cell lysates, tissue homogenates, serum, plasma, urine, as well as other biological fluids when applicable. Run proper controls as necessary. Always run a standard curve with samples. Use 1X Sample Diluent for dilutions and preparations of samples. Sample concentrations must be adjusted to 1-10 mg/mL prior to performing the assay.

Note: For protein samples lower than 1 mg/mL, concentrate by centrifugal filtration or by mixing 1 vol of protein sample and 4 vol of 20% TCA solution, incubating 10 minutes on ice and centrifuging at 10,000 g. Dissolve the protein pellet in a small volume of 1X PBS.

(Optional): High nucleic acid can erroneously contribute to higher estimation of carbonyls. To remove nucleic acid, add streptomycin sulfate or PEI to a final concentration of 1% and 0.5% respectively,

incubate samples 30 minutes at room temperature and remove the nuclei acid precipitates by centrifuging at 6000 g for 10 minutes at 4°C. Test supernatant accordingly.

- Cells or Tissues: Centrifuge cells at 1,500 x g for 5 minutes at 4°C. Remove media and resuspend in 1-2 mL cold 1X Sample Diluent. Homogenize or sonicate cells or tissues on ice. Centrifuge at 10,000 x g for 5 minutes at 4°C. Remove supernatant and store on ice. Store on ice or at -80°C for long-term storage. Prior to use, adjust the total protein concentration to 1-10 mg/mL with 1X Sample Diluent.
- Serum: Collect blood without using an anticoagulant. Allow blood to clot for 30 minutes at room temperature. Centrifuge blood at 2,000 x g for 15 minutes at 4°C. Carefully remove the top yellow serum layer without disturbing the white buffy layer. Store on ice or at -80°C for long-term storage. Prior to use, adjust the total protein concentration to 1-10 mg/mL with 1X Sample Diluent.
- Plasma: Collect blood using an anticoagulant such as citrate or heparin. Centrifuge blood at 1000 x g for 5 minutes at 4°C. Carefully remove the top yellow serum layer without disturbing the white buffy layer. Store on ice or at -80°C for long-term storage. Prior to use, adjust the total protein concentration to 1-10 mg/mL with 1X Sample Diluent.
- Urine: Urine normally has very low protein concentrations and is only recommended if it can be prepared in the 1-10 mg/mL range. Use 1X Sample Diluent for all dilutions.

Preparation of the Fluorophore Standard Curve

1. Use the Fluorophore Standard to prepare a series of standards. Prepare standards just prior to use. Prepare a 25 µM solution of the fluorophore standard by diluting the stock 1:800 with 1X Assay Diluent. Vortex thoroughly. For example, add 2 µL of the 20 mM Fluorophore Standard stock to 1598 µL of 1X Assay Diluent.
2. Prepare a dilution series of fluorophore standards in the concentration range of 0 nM – 1250 nM by diluting the 25 µM fluorophore solution in 1X Assay Diluent (see Table 1).

Standard Tubes	25 µM Fluorophore (µL)	Assay Diluent (µL)	Fluorophore (nM)
1	50	950	1250
2	500 of Tube #1	500	625
3	500 of Tube #2	500	313
4	500 of Tube #3	500	156
5	500 of Tube #4	500	78
6	500 of Tube #5	500	39
7	500 of Tube #6	500	20
8	0	500	0

Table 1. Preparation of Fluorophore Standards.

Assay Protocol

1. Determine the protein concentration of each sample. Adjust the protein concentration to between 1-10 mg/mL with 1X Sample Diluent.
2. Add 50 μ L of sample to a microcentrifuge tube, and follow with a 50 μ L of the freshly prepared 1X Protein Carbonyl Fluorophore solution to each sample tube. Vortex thoroughly.
3. Incubate samples overnight at room temperature and protected from light.
4. Add 400 μ L of 20% TCA solution to each tube. Vortex thoroughly and incubate on ice for 10 minutes.
5. Centrifuge the tubes at 10,000 x g for 10 minutes. Remove and discard the supernatant.
6. Add 1 mL of acetone to each tube and vortex thoroughly in order to break up the pellet. If the pellet does not disintegrate upon vortexing, use a spatula or other small implement to manually break up the pellet.
7. Centrifuge the tubes at 10,000 x g for 10 minutes.
8. Repeat the acetone wash two additional times.
9. After the last wash, remove and discard the supernatant. Leave sample tube lids open, allow the sample pellet to dry out thoroughly for 1 hour.
10. After ensuring that all of the acetone has evaporated, add 50 μ L of Protein Solubilization Solution to each tube. Vortex thoroughly and incubate 10 minutes at room temperature.
11. Add 450 μ L of diluted Assay Diluent to each tube and vortex. Note that samples have been diluted 1:10 from their original concentration.
12. Centrifuge the tubes at 10,000 x g for 10 minutes to remove excess debris.
13. Determine the protein concentration of each sample.
14. Prepare a standard curve as outlined in Table 1.
15. Transfer 100 μ L of each sample or prepared standards to a 96-well black fluorescence microtiter plate.
16. Read the fluorescence with a fluorescence plate reader with 480 nm excitation /530 nm emission filter settings.

Calculation of Results

The following figures demonstrate typical Protein Carbonyl Fluorometric Assay results. Fluorescence measurement was performed on SpectraMax Gemini XS Fluorometer (Molecular Devices) with a 485/538 nm filter set and 530 nm cutoff. One should use the data below for reference only. This data should not be used to interpret actual results.

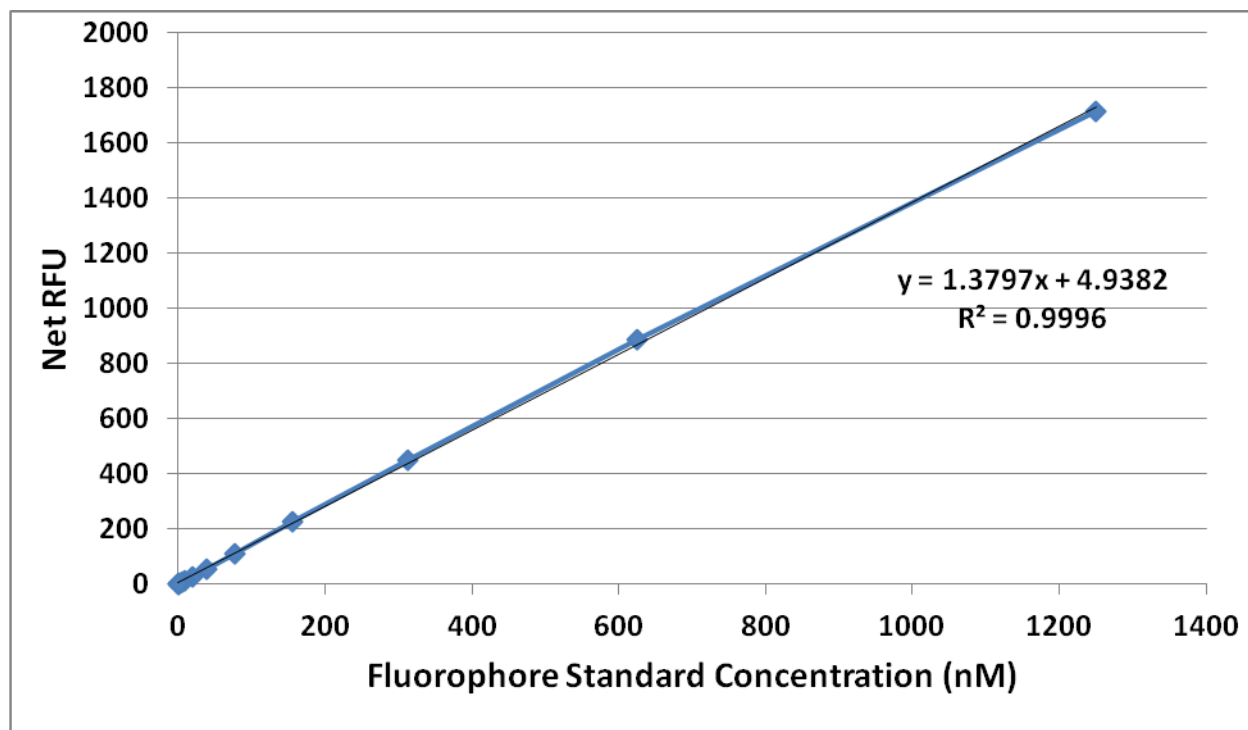


Figure 1. Protein Carbonyl Fluorophore Standard Curve.

Calculations:

1. Determine the average fluorescence for each standard and sample. Subtract the value of 0 nM Fluorophore blank from itself and all other standard and sample values. These are the adjusted RFU values.
2. Graph the adjusted RFU values for the standards as a function of their corresponding concentration from Table 1.
3. Determine the carbonyl concentration of samples using the equation obtained from the linear regression of the standard curve.

Calculation Example:

Chemically oxidized and reduced BSA were prepared at 10, 5, and 1 mg/mL in 1X Sample Diluent. Samples were then derivatized with the Protein Carbonyl Fluorophore and processed according to the Assay Protocol. After TCA precipitation and wash, the protein pellet was dissolved in Protein Solubilization Solution and the protein concentration of the solubilized protein was determined by BCA assay. For example, for the 10 mg/mL Oxidized BSA sample, after 10-fold dilution with Assay Diluent the solubilized protein concentration was 0.63 mg/mL and its adjusted RFU reading is 1015.

$$\text{Protein Carbonyl Concentration (nM)} = [(1015 \text{ RFUs} - 4.9382 \text{ (y intercept)}) / 1.3797 \text{ (slope)}] \times 10 \text{ (Dilution Factor)} = 7321 \text{ nM}$$

$$\text{Protein Carbonyl Content (nmol/mg)} = 7321 \text{ nM} / (0.63 \text{ mg/mL} \times 10 \text{ (Dilution Factor)}) = (7.321 \text{ nmol/mL}) / (6.3 \text{ mg/mL}) = 1.162 \text{ nmol/mg}$$

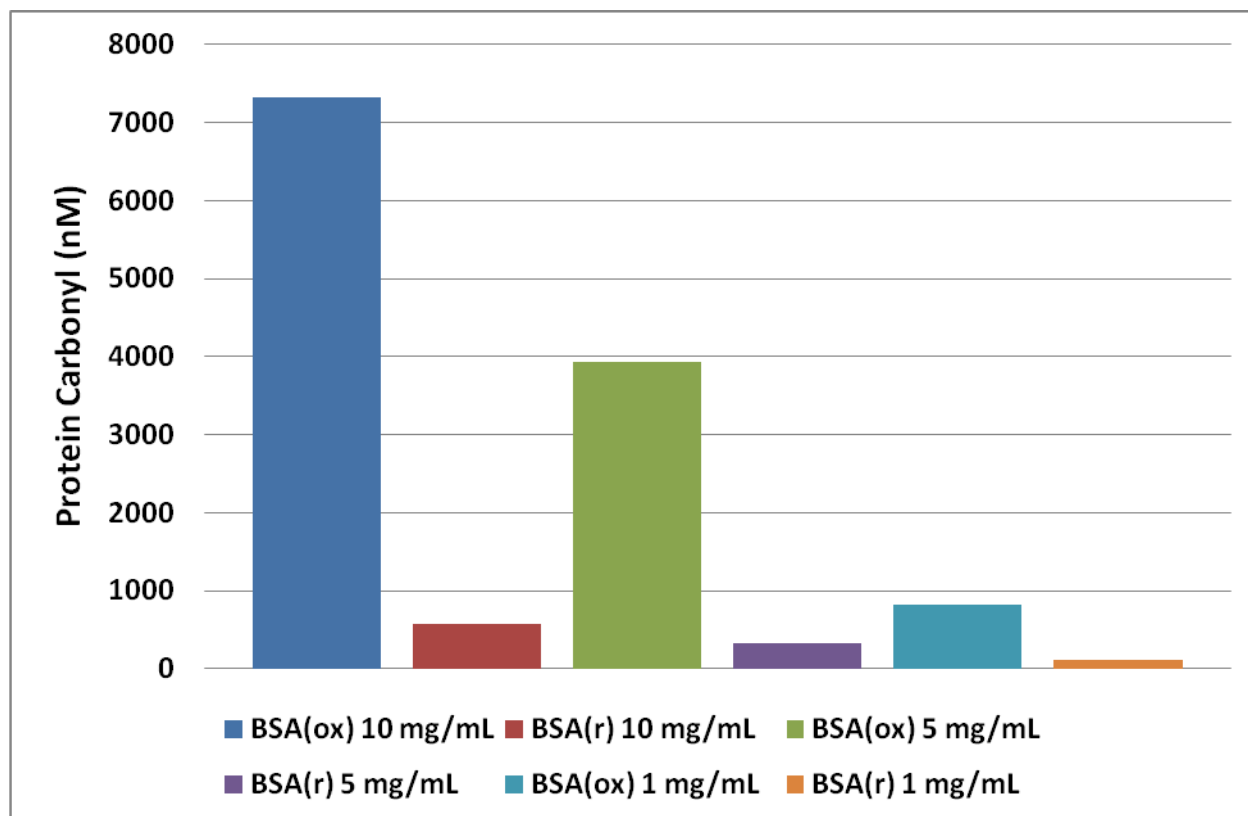


Figure 2. Protein Carbonyl Contents in Oxidized and Reduced BSA.

References

1. Buss, H, Chan, TP, Sluis, KB, Domigan, NM, and Winterbourn, CC. (1985) *Free Radic Biol Med.* **23**: 361-6.
2. Cadenas, E., Boveris, A., Ragan, CI., and Stoppani, AO. (1977) *Archives of Biochemistry & Biophysics* **180**: 248–257.
3. Hammer, M., Schweitzer, D., Richter, S., Konigsdorffer, E. (2005) *Physiol. Meas.* **26**: N9-N12.
4. Reznick, AZ., and Packer, L. (1994) *Methods Enzymol.* **233**: 263-357.
5. Talent, JM., Kong, Y., and Gracy, RW. (1998) *Anal. Biochem.* **263**: 31–38.
6. Wakeyama, H., Takeshige, K., Takayanagi, R., and Minakami, S. (1982) *Biochem J.* **205**: 593–601.

Recent Product Citations

1. Dos Anjos, C. et al. (2023). New Insights into the Bacterial Targets of Antimicrobial Blue Light. *Microbiol Spectr.* **11**(2):e0283322. doi: 10.1128/spectrum.02833-22.
2. Gudiksen, A. et al. (2021). Ameliorating effects of lifelong physical activity on healthy aging and mitochondrial function in human white adipose tissue. *J Gerontol A Biol Sci Med Sci.* doi: 10.1093/gerona/glab356.
3. Nukala, S.B. et al. (2021). Protein network analyses of pulmonary endothelial cells in chronic thromboembolic pulmonary hypertension. *Sci Rep.* **11**(1):5583. doi: 10.1038/s41598-021-85004-z.

4. de Los Santos-Jiménez, J. et al. (2021). Glutaminase isoforms expression switches microRNA levels and oxidative status in glioblastoma cells. *J Biomed Sci.* **28**(1):14. doi: 10.1186/s12929-021-00712-y.
5. Christiansen, L.B. et al. (2021). Atorvastatin impairs liver mitochondrial function in obese Göttingen Minipigs but heart and skeletal muscle are not affected. *Sci Rep.* **11**(1):2167. doi: 10.1038/s41598-021-81846-9.
6. Wake, H. et al. (2020). Histidine-rich glycoprotein possesses anti-oxidant activity through self-oxidation and inhibition of hydroxyl radical production via chelating divalent metal ions in Fenton's reaction. *Free Radic Res.* doi: 10.1080/10715762.2020.1825703.
7. Kim, T. et al. (2020). Cupric ion in combination with hydrogen peroxide and hydroxylamine applied to inactivation of different microorganisms. *J Hazard Mater.* doi: 10.1016/j.jhazmat.2020.123305.
8. Sambon, M. et al. (2020). Dibenzoylethylamine Has Powerful Antioxidant and Anti-Inflammatory Properties in Cultured Cells and in Mouse Models of Stress and Neurodegeneration. *Biomedicines.* **8**(9): E361. doi: 10.3390/biomedicines8090361.
9. Buch, B.T. et al. (2020). Colchicine treatment impairs skeletal muscle mitochondrial function and insulin sensitivity in an age-specific manner. *FASEB J.* doi: 10.1096/fj.201903113RR.
10. Pavlov, D. et al. (2020). Enhanced conditioning of adverse memories in the mouse modified swim test is associated with neuroinflammatory changes - effects that are susceptible to antidepressants. *Neurobiol Learn Mem.* doi: 10.1016/j.nlm.2020.107227.
11. Gorlova, A. et al. (2019). Alteration of oxidative stress markers and behavior of rats in a novel model of depression. *Acta Neurobiol Exp (Wars).* **79**(3):232-237. doi: 10.21307/ane-2019-021.
12. Halling, J.F. et al. (2019). PGC-1 α regulates mitochondrial properties beyond biogenesis with aging and exercise training. *Am J Physiol Endocrinol Metab.* doi: 10.1152/ajpendo.00059.2019.
13. Tiwari, M. K. et al. (2019). Copper ion / H₂O₂ oxidation of Cu/Zn-Superoxide dismutase: Implications for enzymatic activity and antioxidant action. *Redox Biology.* doi: 10.1016/j.redox.2019.101262.
14. Sambon, M. et al. (2019). Thiamine and benfotiamine protect neuroblastoma cells against paraquat and β -amyloid toxicity by a coenzyme-independent mechanism. *Heliyon.* **5**(5), e01710. doi: 10.1016/j.heliyon. 2019.e01710.
15. Gorlova, A. et al. (2019). Thiamine and benfotiamine counteract ultrasound-induced aggression, normalize AMPA receptor expression and plasticity markers, and reduce oxidative stress in mice. *Neuropharmacology.* pii: S0028-3908(19)30058-9. doi: 10.1016/j.neuropharm.2019.02.025.
16. Naparło, K. et al. (2019). Flavanols protect the yeast *Saccharomyces cerevisiae* against heating and freezing/thawing injury. *J Appl Microbiol.* **126**(3):872-880. doi: 10.1111/jam.14170.
17. Frontiñán-Rubio, J. et al. (2018). Sex-dependent co-occurrence of hypoxia and β -amyloid plaques in hippocampus and entorhinal cortex is reversed by long-term treatment with ubiquinol and ascorbic acid in the 3 \times Tg-AD mouse model of Alzheimer's disease. *Mol Cell Neurosci.* **92**:67-81. doi: 10.1016/j.mcn.2018.06.005.
18. Adamczyk-Sowa, M. et al. (2017). Oxidative Modification of Blood Serum Proteins in Multiple Sclerosis after Interferon Beta and Melatonin Treatment. *Oxid Med Cell Longev.* **2017**:7905148. doi: 10.1155/2017/7905148.
19. Yeh, A. et al. (2017). Effect of contaminants of emerging concern on liver mitochondrial function in Chinook salmon. *Aquat. Toxicol.* **190**:21-31.

20. Tangtrongsup, S. and Kisiday, J.D. (2017). Modulating the oxidative environment during mesenchymal stem cells chondrogenesis with serum increases collagen accumulation in agarose culture. *J. Orthop. Res.* doi:10.1002/jor.23618.
21. Adamczyk-Sowa, M. et al. (2017). Oxidative modifications of blood serum proteins in myasthenia gravis. *Journal of Neuroimmunology.* **305**:145-153.
22. Rey, B. et al. (2016). Hormetic response triggers multifaceted anti-oxidant strategies in immature king penguins (*Aptenodytes patagonicus*). *Free Radic Biol Med.* doi: 10.1016/j.freeradbiomed.2016.07.015.
23. Li-Byarlay, H. et al. (2016). Honey bee (*Apis mellifera*) drones survive oxidative stress due to increased tolerance instead of avoidance or repair of oxidative damage. *Exp Gerontol.* doi: 10.1016/j.exger.2016.07.003.
24. Maslanka, R. et al. (2016). The effect of berry juices on the level of oxidative stress in yeast cells exposed to acrylamide. *J Food Biochem.* doi: 10.1111/jfbc.12260.
25. Lauritzen, K. H. et al. (2015). Impaired dynamics and function of mitochondria caused by mtDNA toxicity leads to heart failure. *Am J Physiol Heart Circ Physiol.* **309**:H434-H449.
26. Sadowska-Bartosz, I. & Bartosz, G. (2015). Ascorbic acid and protein glycation in vitro. *Chem Biol Interact.* doi: 10.1016/j.cbi.2015.07.006.
27. Zabala, V. et al. (2015). Potential contributions of the tobacco nicotine-derived nitrosamine ketone (NNK) in the pathogenesis of steatohepatitis in a chronic plus binge rat model of alcoholic liver disease. *Alcohol Alcohol.* doi: 10.1093/alcalc/agu083.
28. Rey, B. et al. (2014). Thyroid status affects membranes susceptibility to free radicals and oxidative balance in skeletal muscle of Muscovy ducklings (*Cairina moschata*). *J Exp Zool A Ecol Genet Physiol.* **321**:415-421.
29. Tong, M. et al. (2014). Therapeutic reversal of chronic alcohol-related steatohepatitis with the ceramide inhibitor myriocin. *Int J Exp Pathol.* **95**:49-63.

Warranty

These products are warranted to perform as described in their labeling and in Cell Biolabs literature when used in accordance with their instructions. THERE ARE NO WARRANTIES THAT EXTEND BEYOND THIS EXPRESSED WARRANTY AND CELL BIOLABS DISCLAIMS ANY IMPLIED WARRANTY OF MERCHANTABILITY OR WARRANTY OF FITNESS FOR PARTICULAR PURPOSE. CELL BIOLABS' sole obligation and purchaser's exclusive remedy for breach of this warranty shall be, at the option of CELL BIOLABS, to repair or replace the products. In no event shall CELL BIOLABS be liable for any proximate, incidental or consequential damages in connection with the products.

Contact Information

Cell Biolabs, Inc.
 7758 Arjons Drive
 San Diego, CA 92126
 Worldwide: +1 858-271-6500
 USA Toll-Free: 1-888-CBL-0505
 E-mail: tech@cellbiolabs.com
www.cellbiolabs.com

©2011-2023: Cell Biolabs, Inc. - All rights reserved. No part of these works may be reproduced in any form without permissions in writing.