

Complete Classic Medium Kit With Serum and CultureBoost™

Cat # NB-11-0046

Description

This medium is formulated with 10% FBS serum. This becomes a complete medium once activated with the included CultureBoost™ growth factor supplement. Basal medium contains no added hormones and contains no antibiotics. This medium offers consistent and stable environment to optimize the thawing and passaging of primary cells. The serum-based formulations:

- Are based on a modified and optimized basal DMEM/F12 (1:1)
- Contain serum at 10% by volume
- Do not contain phenol red
- Packaged in 500mL bottles

Bac-Off® (Cat # NB-11-0074) antibiotic solution is recommended for all Neo Biotech media. Medium contains no phenol red, which interferes with cell physiology.

Neo Biotech media and reagents are sterile, made with WFI and all components are cGMP and ISO Compliant.

Components

Kit consists of 500mL Medium, 5mL CultureBoost™ (Cat # NB-11-0066, containing animal derived growth factors) and 10mL Attachment Factor™ (Cat # NB-11-0088, extracellular matrix reagent for coating flasks).

Storage & Shipping

Immediately refrigerate (2 - 8°C) the medium, CultureBoost™, and AttachmentFactor™. If the entire 500mL unit of media will not be used within 30 days, supplement the medium with CultureBoost™, then aliquot and freeze in smaller units, which must be used within 30 days. Once the unit is supplemented, or any component of the medium kit is opened, the shelf life is 30 days at refrigerated (2 - 8°C) temperatures. Store Attachment Factor™ at 2 - 8°C between uses.

All medium and reagents shipped at ambient temperature.

Product Use:

NB-11-0046 is for laboratory research only. It is not approved for human or animal use, or for application in *in vitro* diagnostic procedures.

Instructions for Use

Spray the outside of all tools/instruments (bottles, tubes, racks, etc.) with ethanol solution before bringing them under the hood. Make sure all work surfaces are disinfected. Spray the outside of your gloves prior to work under the hood.

- Place under hood 500mL media and 5mL CultureBoost™
- Remove shrink wrap and place cap sides down under hood onto ethanol sprayed gauze.
- Pipette or pour 5mL CultureBoost™ into 500mL media.
- Replace cap on media and refrigerate or use in cell culture.

Media is now complete.

***For reference only
For Research Use Only. Not for Diagnostic or Therapeutic Use.***