

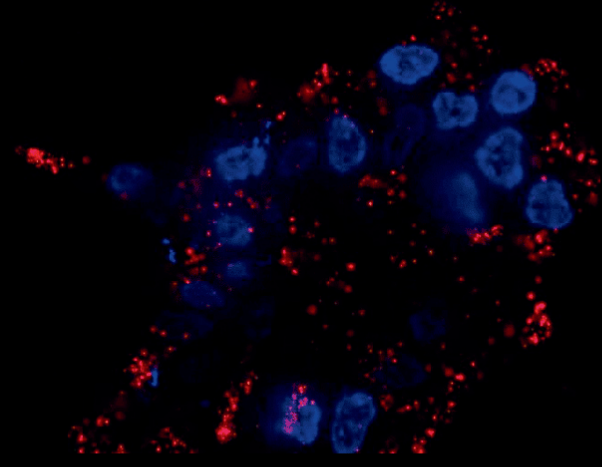
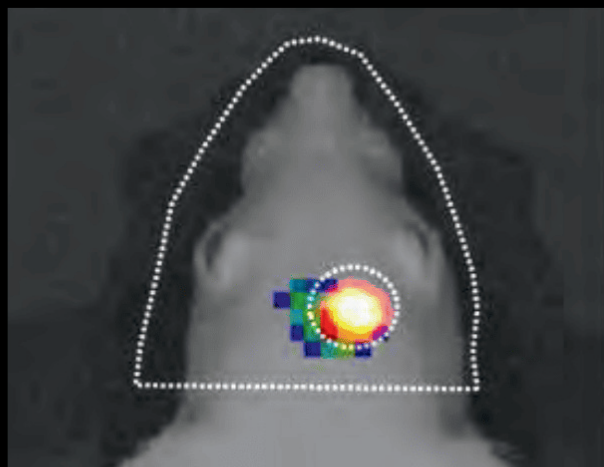
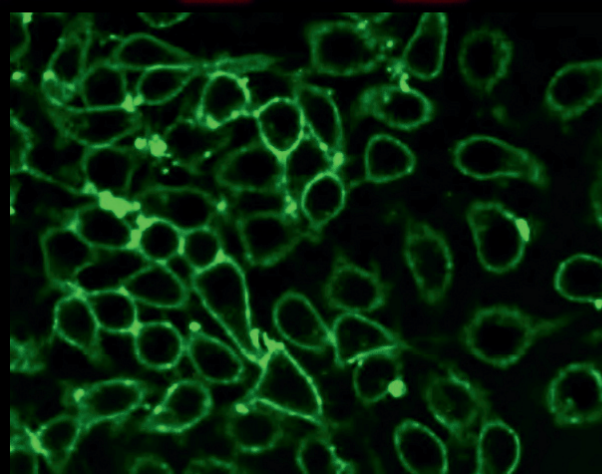


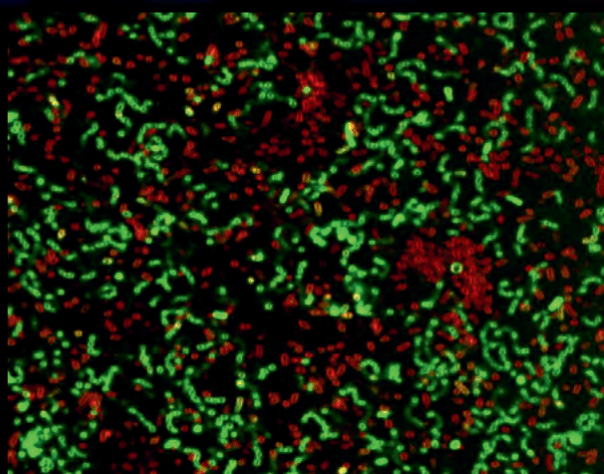

Empowering Life Science with Precision

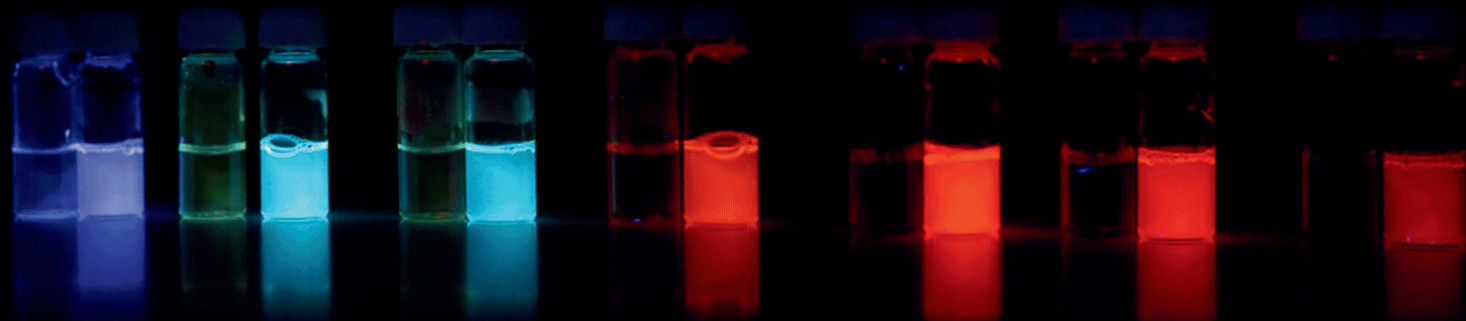

Optical Probes for Mammalian Cells, Extracellular Vesicles & More

Technical Bulletin

Distributed by

cliniSciences



JOTDYES IN A SNAPSHOT

02

Innovative dye technology

MAMMALIAN CELLS

03

Membrane dye| Cell division tracking| RBC

LIPOSOMES

05

Single particle analysis

EV RESEARCH

08




Detection| Isolation| Tracking

JotDyes in a Snapshot



Tech at a Glance

- Workflow compatible
- Accurate
- Variety of Colors (UV to NIR)
- Water Soluble
- Innovative Chemistry
- "Light up" Mechanism

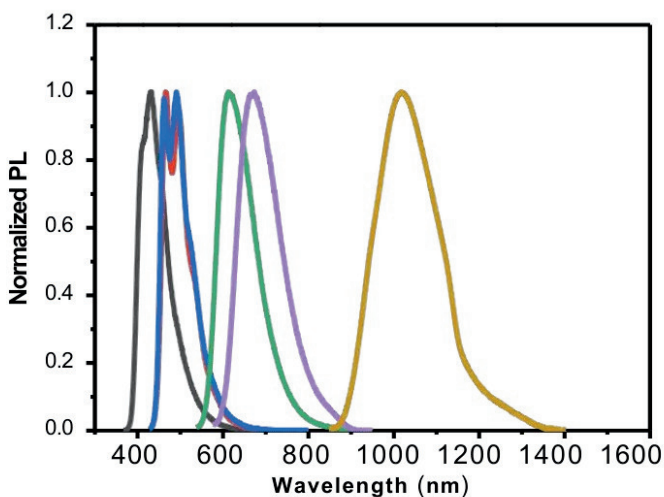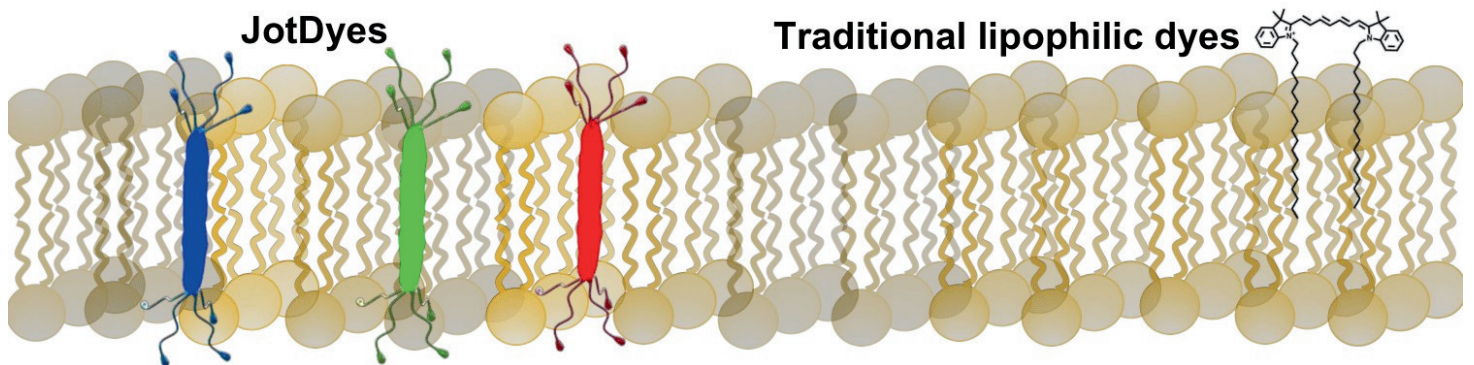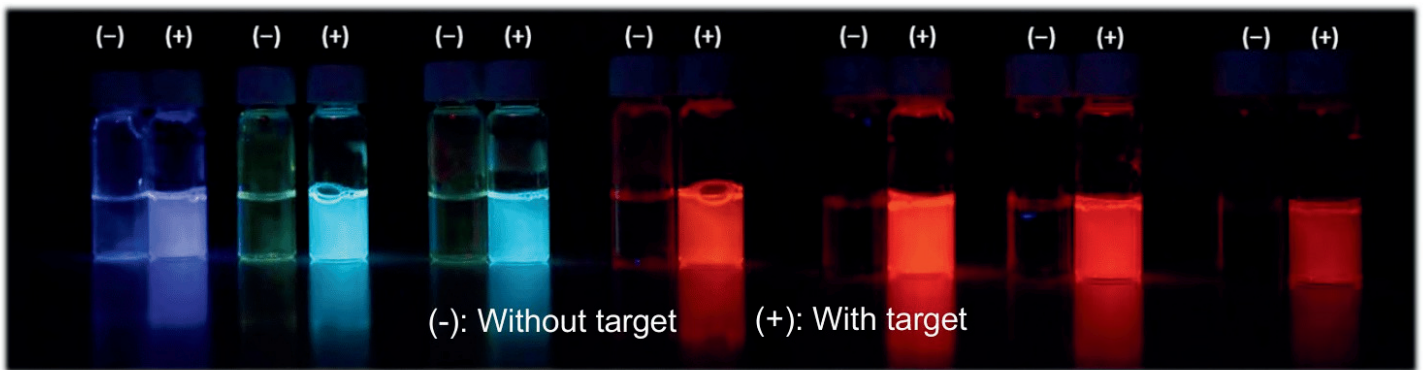



With a variety of options to fit your needs, Jotbody provides a novel class of revolutionary dyes that will **accelerate your research**



Jotbody Membrane Labeling Dye

Confocal images of HepG2 cells that were co-stained with Jotbody Membrane Labeling Dye and a commercially available membrane dye, FM 4-64.

(a) JotDye channel (Ex405), (b) FM 4-64 channel (Ex488), (c) brightfield channel, (d) merged channel. The co-localization of Jotbody Membrane Dye and FM4-64 demonstrates the membrane specificity of JotDye.

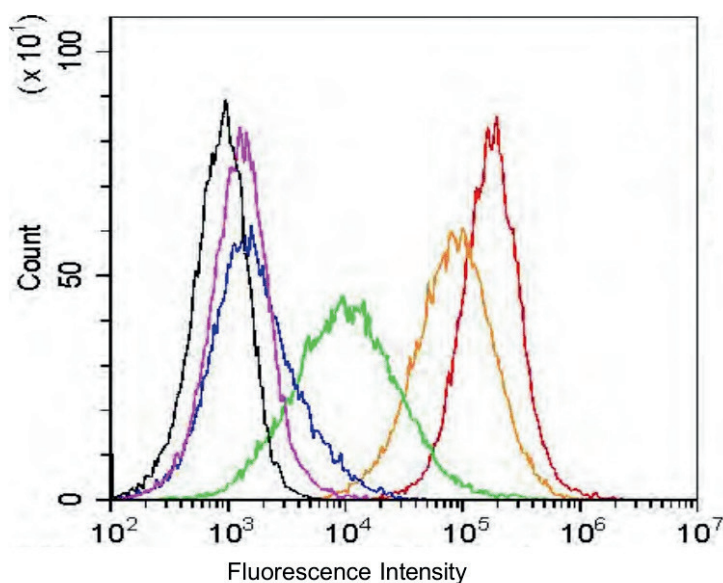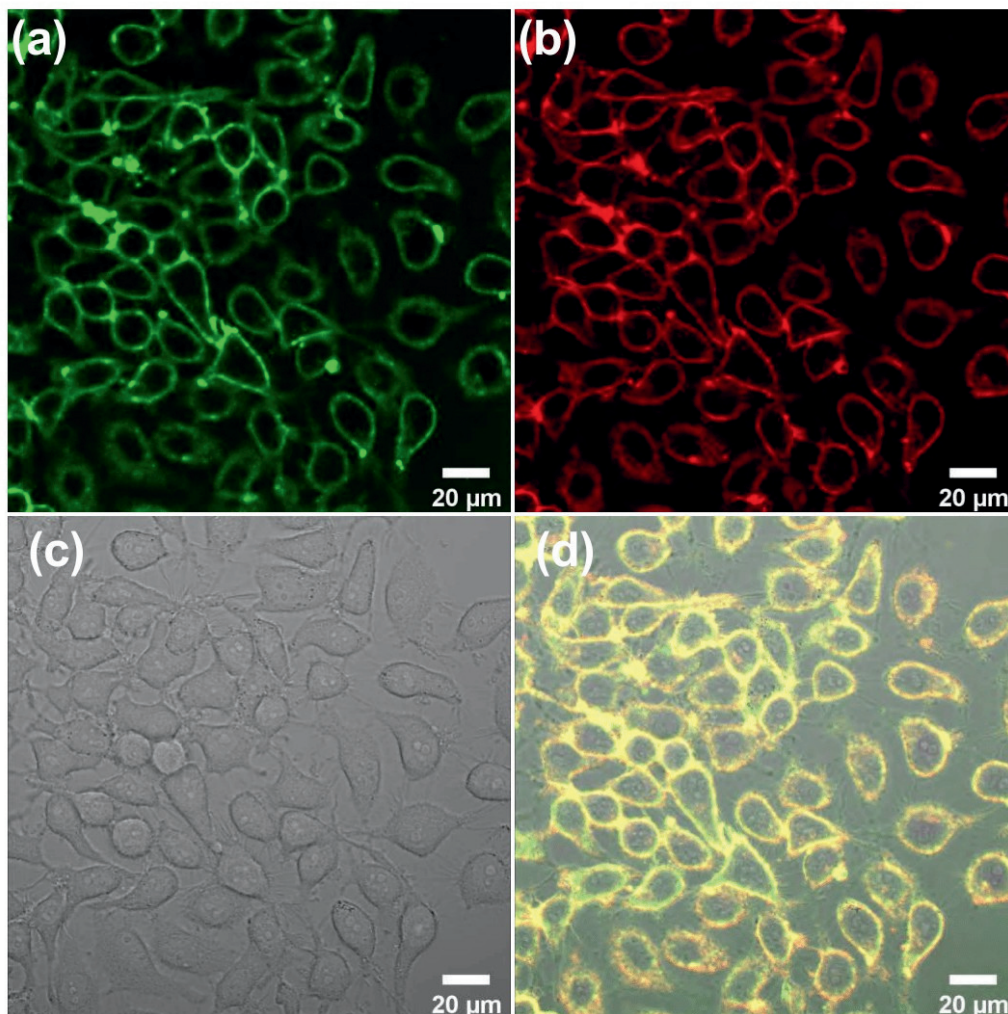


JotDye for Cell Division Tracking

A549 mammalian tumor cells labeled by Jotbody Cell Division Tracking dye. JotDye-labeled cells (red) have a 3-log fold increase in fluorescence intensity compared to the unstained cells (black) were maintained over 4 passages through serial passaging before analysis on CytoFLEX (Beckman Coulter). Orange: 1st passage, Green: 2nd passage, Blue: 3rd passage. Magenta: 4th passage



JotDye for Labeling Red Blood Cells

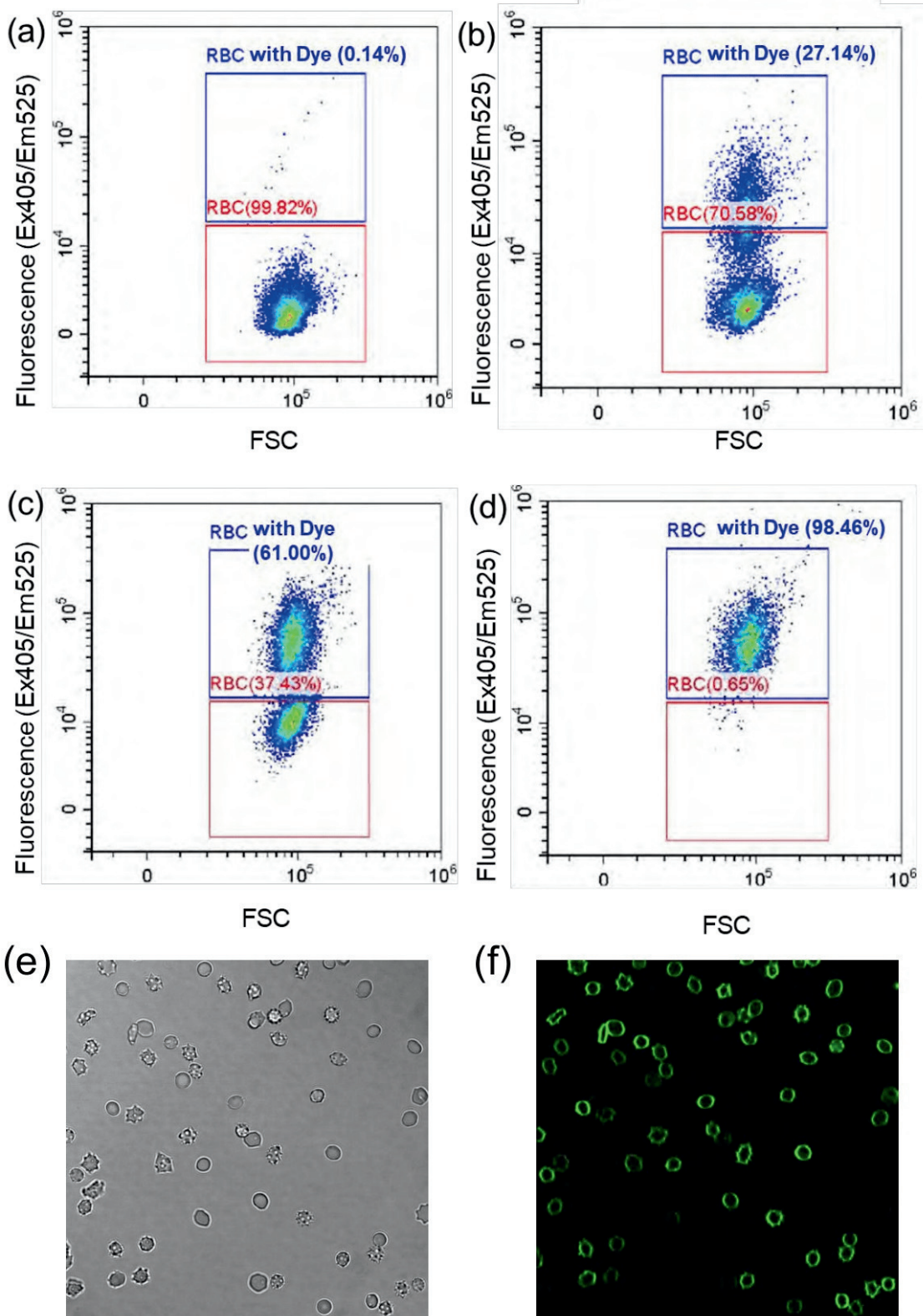


Red blood cells. Flow cytometry analysis of samples that have a mixture of unstained red blood cells (RBC) and RBC stained with JotDye, in the ratio of (a) 0:5, (b) 2:3, (c) 3:2 and (d) 5:0. The JotDye for RBC remains stable in the cell after labelling and does not cross-over into samples that are not intended to be labelled with the dye. (e) Bright field and (f) fluorescent confocal micrograph of RBCs stain with JotDye.

Distributed by

clinisciences




Jotbody Membrane Dyes: Colocalization with traditional lipophilic dyes

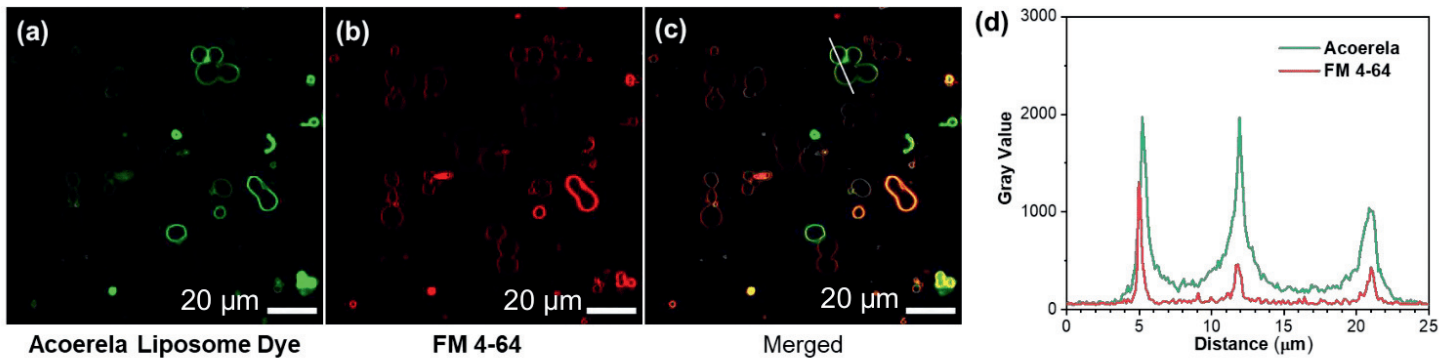

Localization of Jotbody Liposome Dye in the lipid bilayer. Large multilamellar vesicles, comprising 1- palmitoyl-2-oleoyl-glycero-3-phosphocholine (POPC) and 1-palmitoyl-2-oleoyl-sn-glycero-3-phospho-1'- glycerol (POPG) (8:2 ratio), were stained simultaneously with 15 μM of Jotbody Liposome Dye and FM4-64. Fluorescence channels for (a) JotDye, (b) FM4-64 and (c) both channels merged. (d) Fluorescence trace for the white line in (c) demonstrates the colocalization of JotDye and FM4-64.

Pain points with Current Lipophilic Dyes

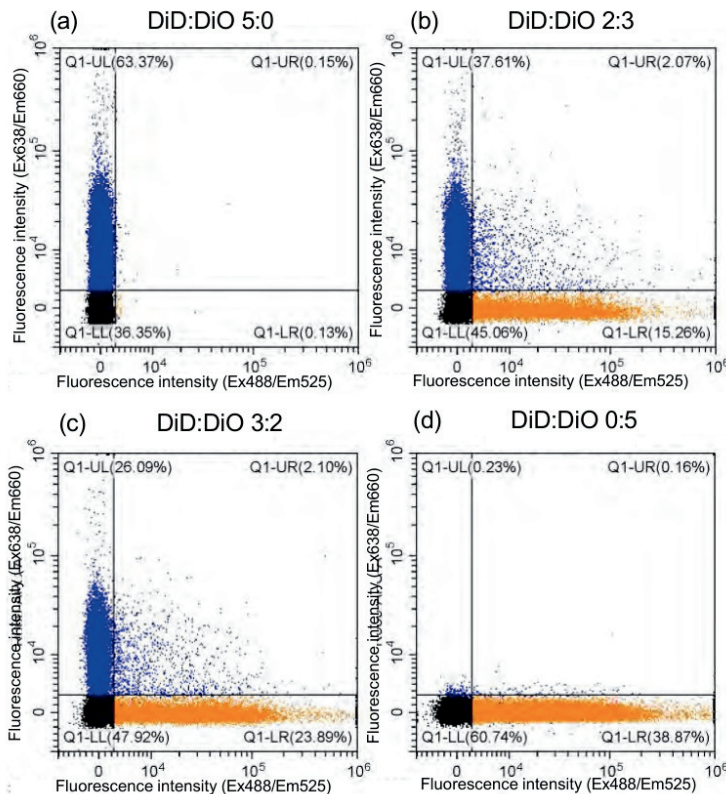


Instability of liposome labeling using the commercially available carbocyanine dyes. Mixtures of labeled 130 nm liposomes in the ratio of DiD:DiO (a) 0:5, (b) 2:3, (c) 3:2 and (d) 5:0, were analyzed on CytoFLEX. (Beckman Coulter) after 1 h incubation. The presence of dual-positive events in Q1UR suggests that the dye labels were not stable and can transfer from one population to another.

Labeling using lipophilic dyes from the carbocyanine family tend not to be stable due to the "leakage" of dye from their targets. These free dyes end up labeling other membrane-based targets that can lead to misleading conclusions.





Single Particle Analysis using JotDyes Stable labeling: No inter-particle dye exchange

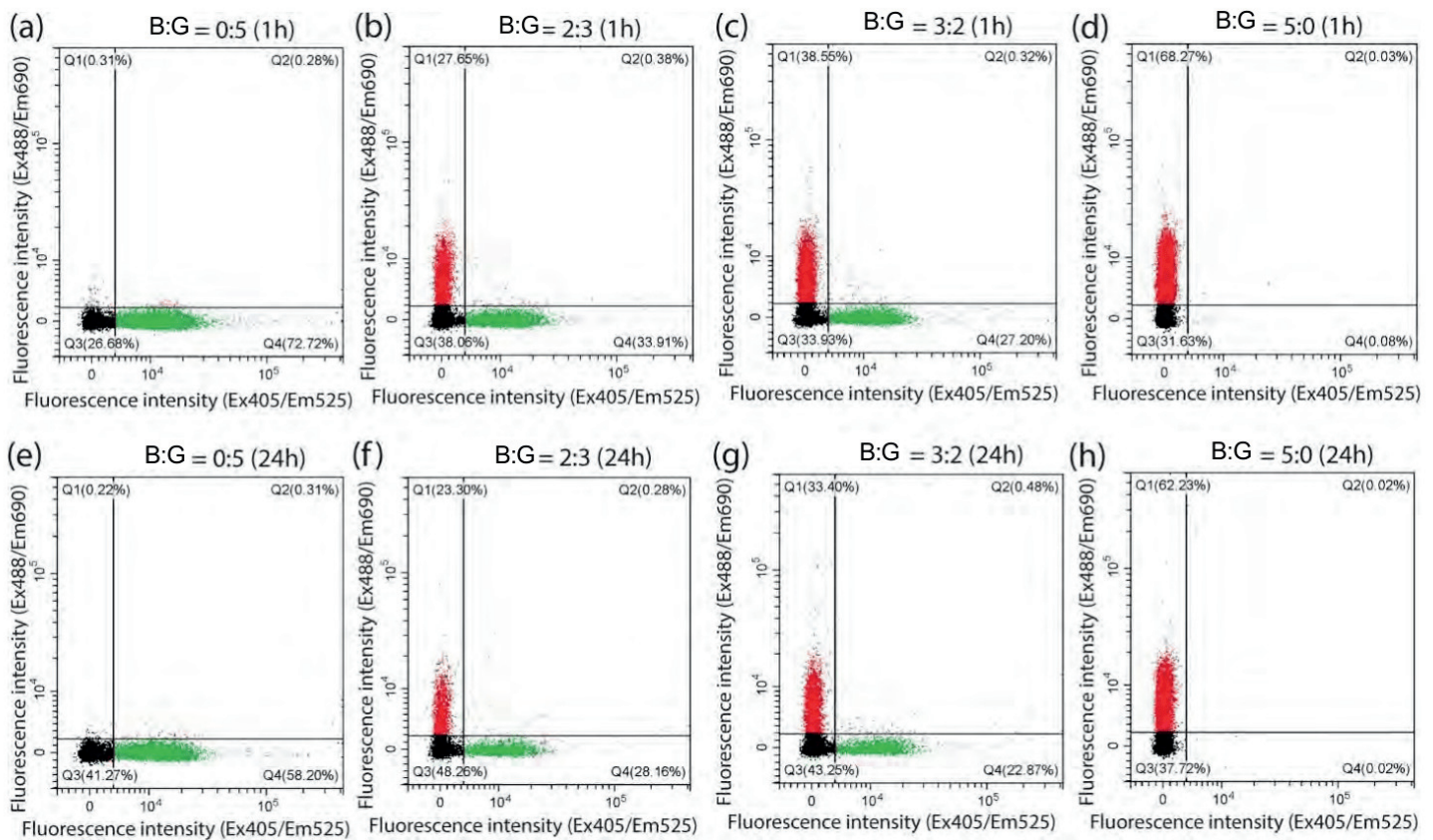

Single particle analysis of 130 nm liposomes labeled with Membrane-Stable Jotbody Blue Dye (B) or Green Dye (G). Mixtures of labeled liposomes in the ratio of Jotbody B:R (a,e) 0:5, (b,f) 2:3, (c,g) 3:2 and (d,h) 5:0, were analyzed on CytoFLEX. (Beckman Coulter) after 1 h (a-d) and 24 h (e-h) incubation. The stability of labeling with JotDyes allowed for a consistent analysis of the labeled liposome mixture even after 24 h, where no dye exchanges were observed between the two liposome populations. The advantage of such stability in dye labeling allows users to unambiguously make conclusions on their observations of the fluorescence data.



No Micelle Formation of JotDyes

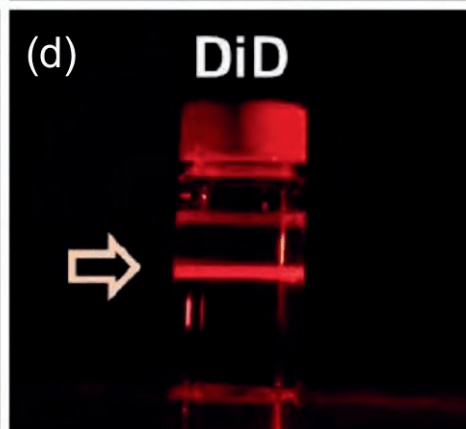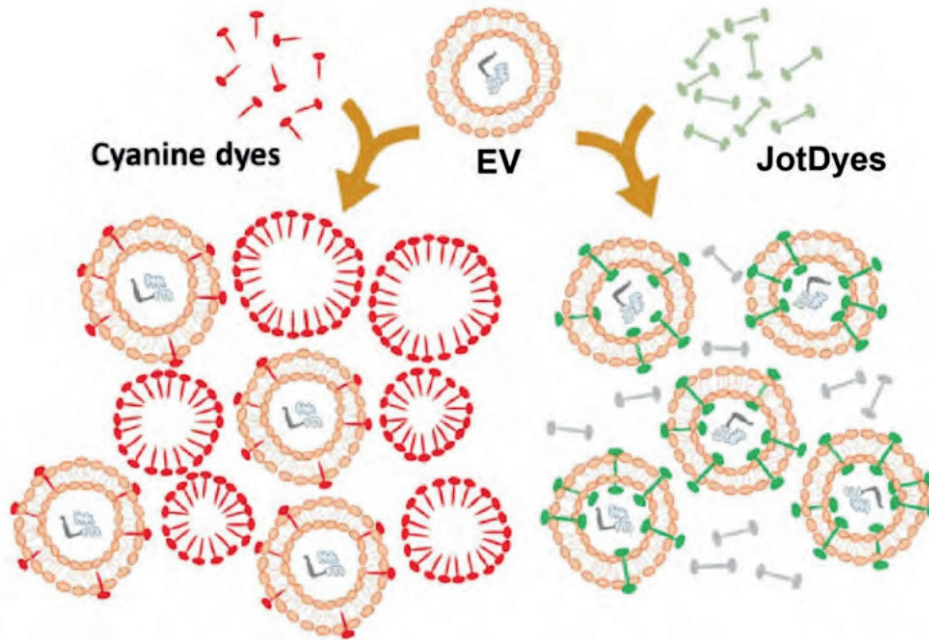


Tyndall effect of nanoparticles. JotDyes are highly water soluble and do not form micelles or nanoparticles when reconstituted in aqueous buffer (a-b), unlike that of the family of PKH26, DiR and DiD dyes (c-d), in which their nanoparticles will give rise to false positives in flow cytometry analysis.

JotDyes promise a **low rate of false positives** for the accurate and reliable detection of EVs.



Using JotDyes for Single Particle Analysis by Flow Cytometry

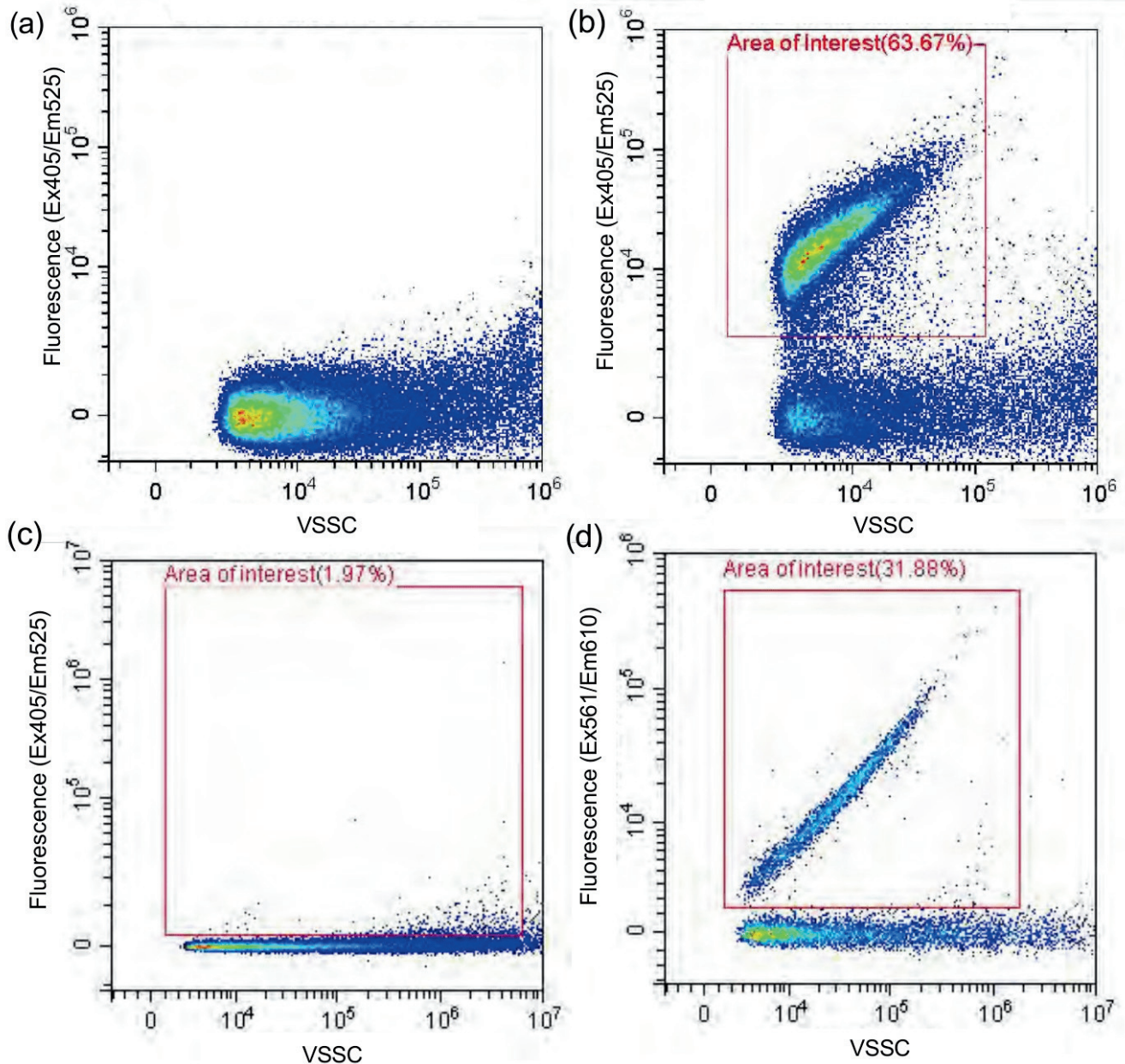


Flow cytometry analysis of EVs labeled with JotDye. Dot plot of (a) unstained EVs, (b) EVs labeled with 1 μ M JotDye, (c) 1 μ M JotDye only and (e) 1 μ M of commercially available PKH26 dye only i.e., in the absence of EVs.

Exosomes are extracellular vesicles that serve as biomarkers for disease detection and have also been utilized as delivery vehicles for therapeutics.





Using JotDyes to Normalize Data for Population Analysis Using Flow Cytometry

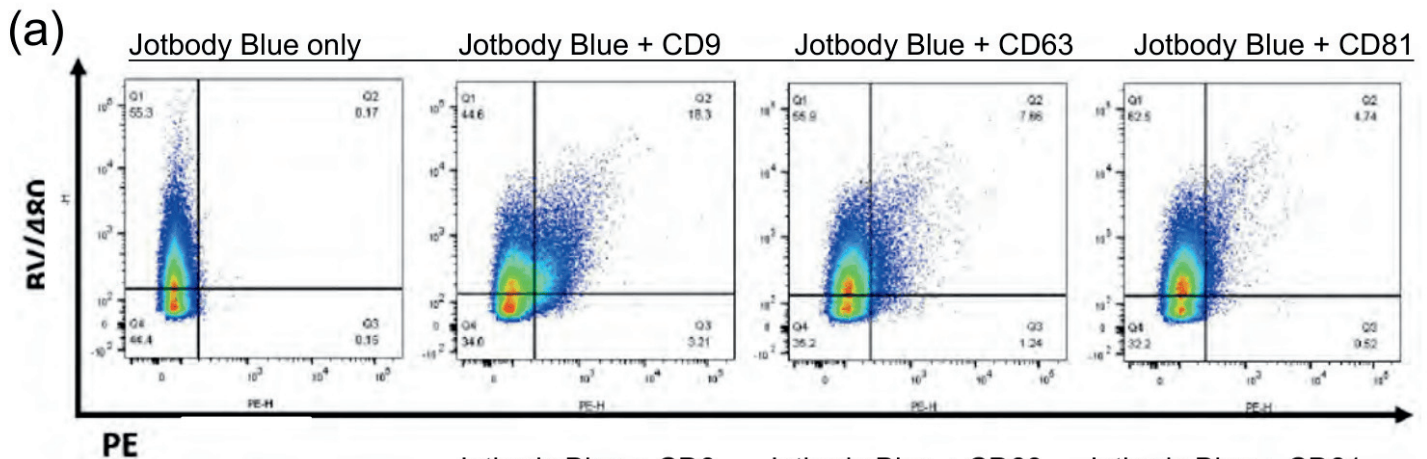

(b)


Gating the Jotbody Blue positive events to normalize the 100% EV population

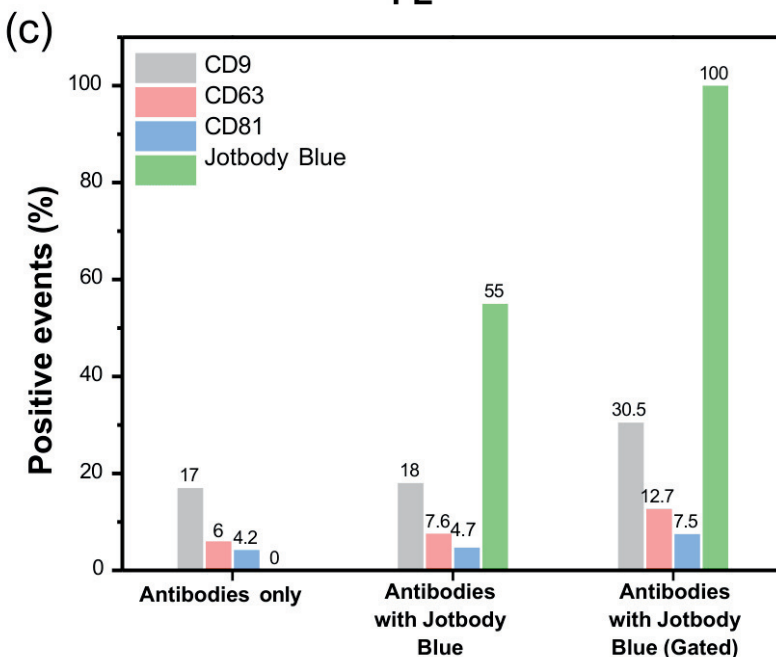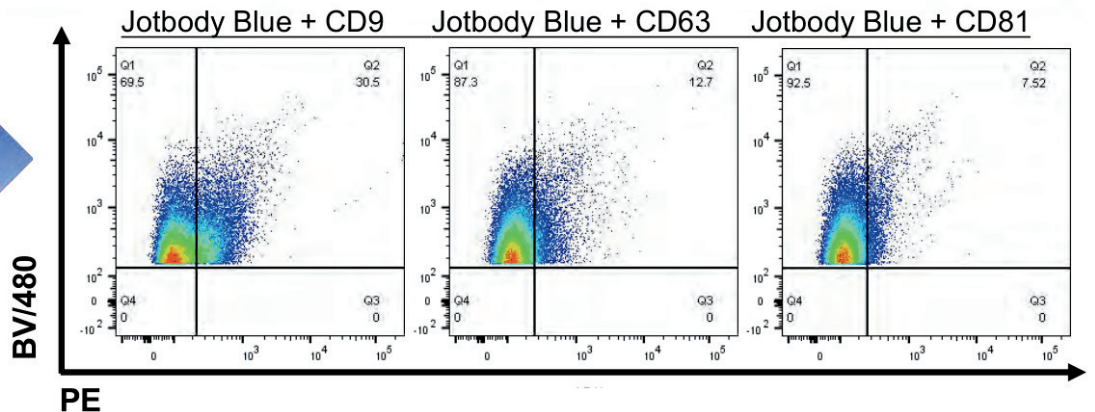

Population analysis of EVs using JotDyes. Flow cytometry measurements were carried out on the A1 BDflow cytometer (Becton Dickinson). (a) Ungated dot plot with x-axis for PE- conjugated antibody channel (PE) and y-axis for Jotbody Blue Dye channel (BV480). From left to right: EVs labeled with Jotbody Blue Dye only, EVs labeled with Jotbody Blue Dye and CD9-PE, with CD63-PE and CD81-PE. (b) The plots in (a) were then gated to only show the Jotbody Blue Dye positive events (i.e., 100% EV population) to reveal the true percentage population of the CD- positive EVs. (c) Plot of dye positive events (%) for EVs that were only stained with antibodies or co-stained with Jotbody Blue Dye, and the effect of gating on the Jotbody positive events.




JotDye-Labeled EV Detection on Nano Flow Cytometry

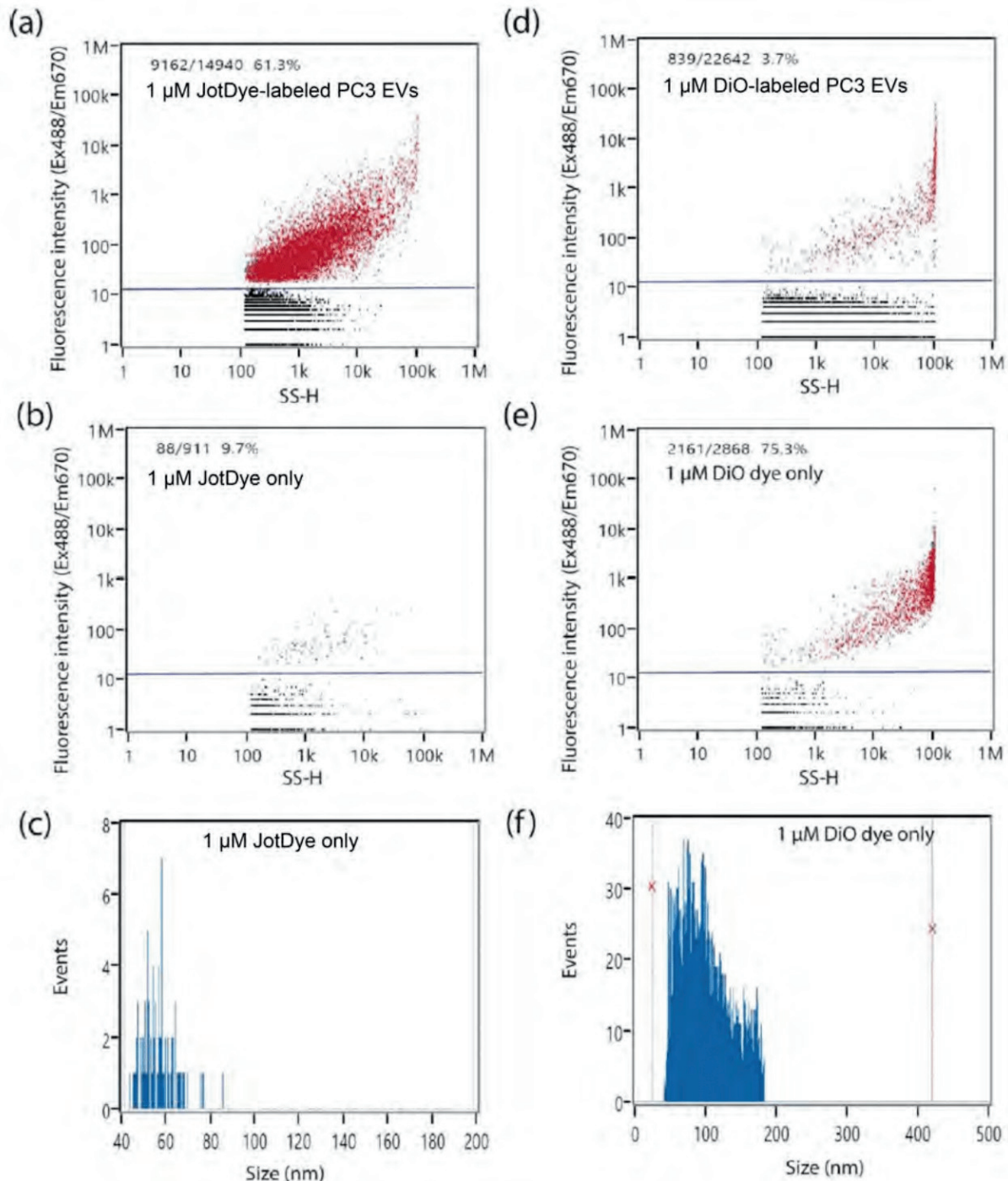

Nano flow cytometry. PC3 EVs were labeled with (a) 1 μ M JotDye and analyzed on the NanoAnalyzer (NanoFCM). A low rate of false positives with the JotDye in buffer only (b) with negligible particle counts detected (c). In contrast, 130 nm-size nanoparticles were observed in a solution of DiO dye only, which did not stain the EVs efficiently (d-f).




Monitoring Uptake of JotDye-Labeled EVs with Fluorescent Microscopy

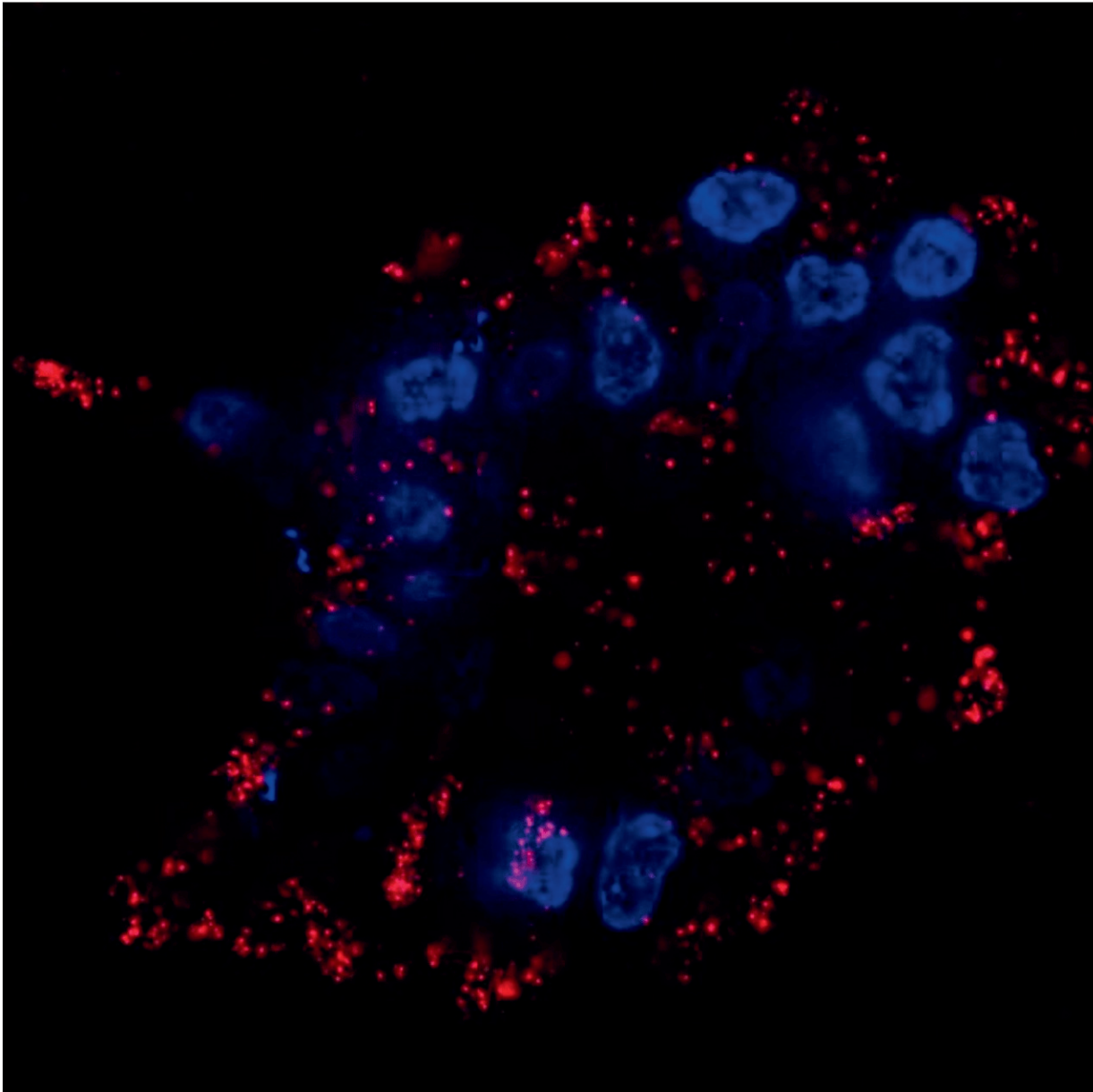

Uptake of JotDye-Labeled EVs in cell culture. Fluorescence micrograph of HEK-293T cells treated with 20 ug of PC-3 EVs stained with JotDye-600 (red) after 24 hours of incubation. Nuclei stained with SYTO Far red (blue). Image taken on Leica THUNDER imager fluorescence microscope and processed using the built-in THUNDER computational clearing method and ImageJ.




Monitoring Uptake of JotDye-Labeled EV with Imaging Flow Cytometry

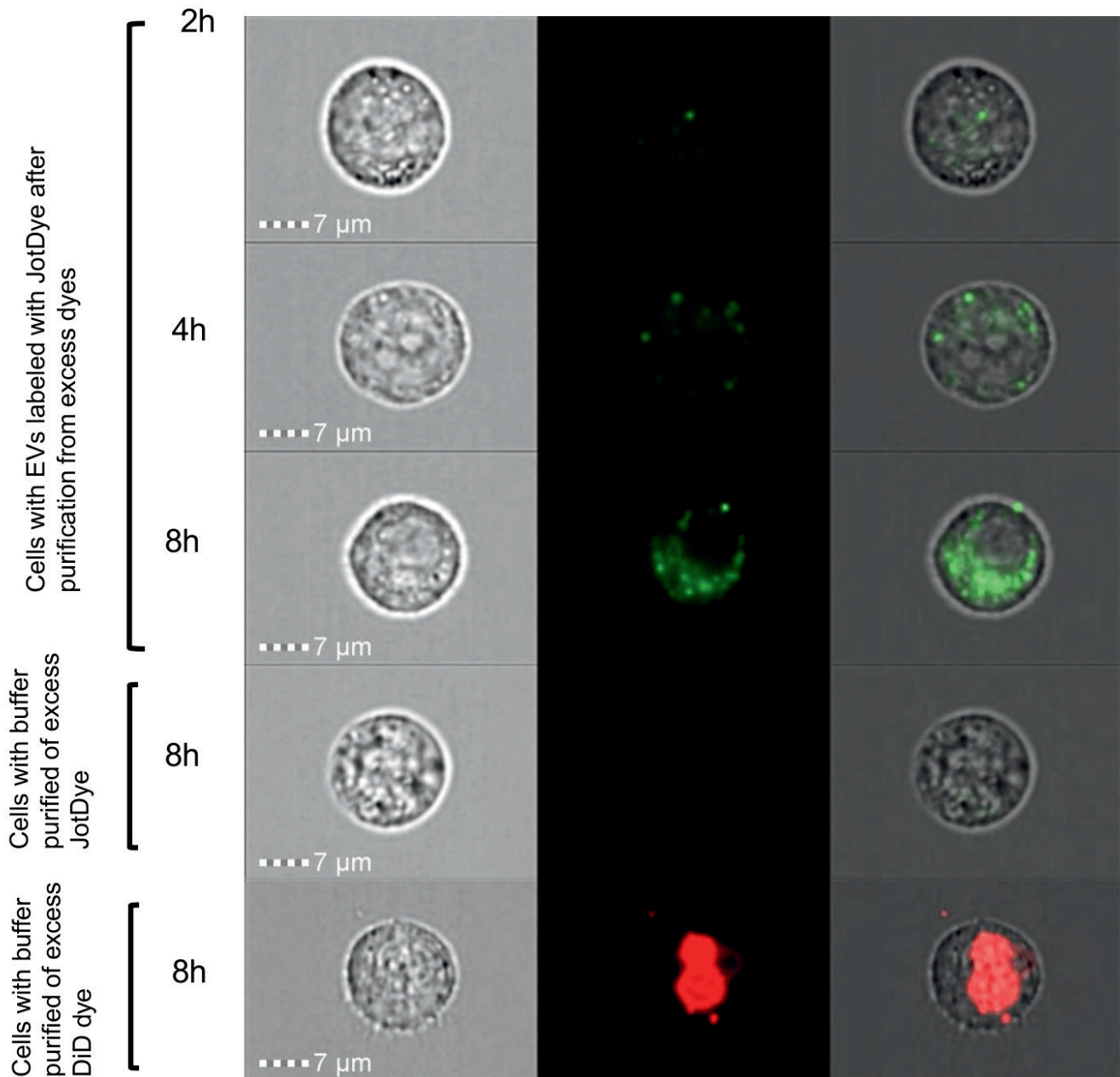

Imaging flow cytometry. PC3 EVs were labeled with 2 μM JotDye and purified using the Amicon 100KDa filter before they were added to A549 cells and incubated. The cells were analyzed on the Imagestream (Amnis) after (a) 2 h, (b) 4 h and (c) 8 h of incubation. The JotDye-labeled EVs were observed as spots (green). In the absence of EVs, no fluorescence signal was observed in the cells because all the free JotDye, which does not form micelles, were removed from the purification step (d). By contrast, micelles of DiD dye were not purified and thus there was still fluorescence signals being observed from the cells, suggesting the uptake of the DiD micelles which can be mistaken for DiD-labelled EVs instead.




JotDye for Endogenously Stained EVs

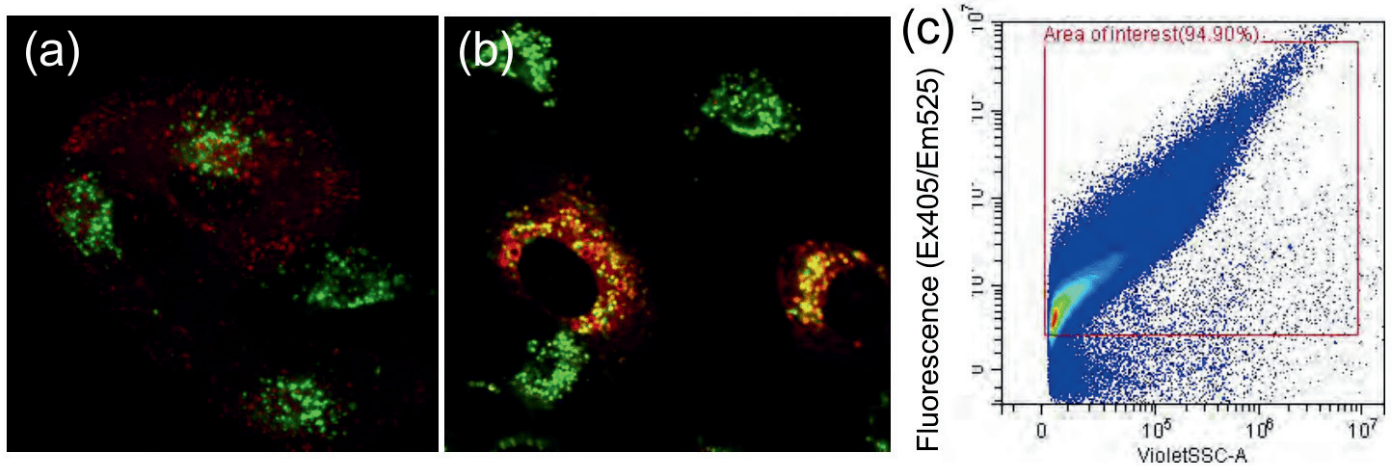

Endogenously stained EVs with JotDye. A549 cells were incubated with 2 μM JotDye for 2 days before the cell culture was ultra-centrifuged to obtain the pre-stained EVs. Confocal imaging studies reveal that the internalized JotDye (green in both (a) and (b)) did not co-localize with the (a) early endosome marker (red) but co-localized with the (b) late endosome marker (red) instead. Late endosomes are the intracellular origins of EV production. (c) Dot plot of stained EVs with JotDye.

TIRF Imaging of JotDye-Labeled EVs

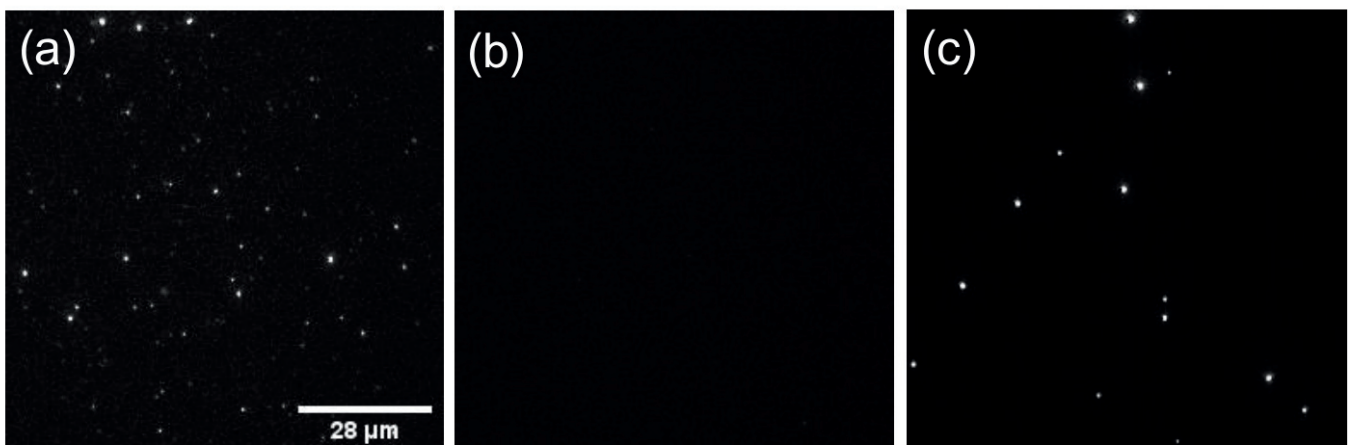

Total internal reflection fluorescence (TIRF) imaging of EVs. PC3 EVs labeled with Jotbody EV Green Dye were immobilized on a substrate and imaged using TIRF microscopy; the diffraction limit is approximately 100 nm. (b) Jotbody EV Green Dye itself does not form any nanoparticles that are of the same size as the EVs. Hence, no fluorescence was observed. On the contrary, the commercially available PKH26 dye forms particles that are large enough to be imaged similarly like the EVs by TIRF (c).

How to navigate through our website?

Homepage: When you first visit our website, you'll land on the homepage. This is where you can find an overview of our services, products, or any important announcements. From here, you can navigate to other sections of the website.

Consult our catalog and services

Choose your Country

Log in to your account

Choose your language

View Cart

Search for a product by description, reference, CAS, Number, target...

upcyte® cells
Expanded primary cells

Categories

- Cell-based assay kits
- Culture flasks
- Culture supports for microscopy
- Detection kits for IHC
- FPLC Columns
- Histological special stains
- Immunoassays
- Lectins
- Next Generation Sequencing (NGS)
- Plant tissue culture
- Radiochemicals
- More categories...

All our products >

Search Bar: Look for specific Products and equipment

Search for a product (description, reference, CAS number, target ...)

Search result : 39126007 product found

Search by selecting either RUO or IVD products

Refine your search by choosing from the following options

Apply Filter

Refine your search :

- RUO
- CE / IVD

Conjugate (100) +

Reactivity species (100) +

Class (66) +

Host species (100) +

Cell name (2) +

Application (100) +

Clone (100) +

APPLY FILTERS

REINITIALIZE

SHARE THIS SEARCH

| Cat# | Description | Cond. |
|-----------------|--|-----------|
| NB-45-00042-100 | Super NI-NTA Agarose Resin | 100ml |
| NB-45-00042-25 | Super NI-NTA Agarose Resin | 25ml |
| NB-45-00058-4 | Proteus 1 -step Batch Mini Spin Column Pack | 40pc |
| NB-12-6001-3 | NeoLine pipette 2-20 µl | 1unit |
| NB-12-0023C | Mini Centrifuge N500C @10,000rpm (Including 6x1.5/2.0ml angle rotor) | 1pcs |
| NB-03-0160 | Proteinase K (Powder) | 100mg |
| NB-60-0001 | NeoPrep mini | 50columns |
| NB-12-8001-19 | Combs for NeoPRO4 mini (1.5mm, 15 wells) | 5pieces |
| NB-12-8001-20 | Spacer glasses flat for NeoPRO4 mini (0.75mm, 100*83mm) | 5pieces |
| NB-12-8001-04 | Short glasses flat for NeoPRO4 mini (1.0mm, 100*73mm) | 10pieces |

“Remember, if you ever get lost or have questions, don’t hesitate to reach out to our customer support team. They’re here to help you address any concerns you may have.”

clinisciences

Austria

Company: CliniSciences GmbH
Address: Sternwartestrasse 76, A-1180
Wien - Austria
Telephone: +43 720 115 580
Fax: +43 720 115 577
Email: oesterreich@clinisciences.com
Web: <https://www.clinisciences.com>



Belgium

Company: CliniSciences S.R.L
Address: Avenue Stalingrad 52, 1000
Brussels - Belgium
Telephone: +32 2 31 50 800
Fax: +32 2 31 50 801
Email: belgium@clinisciences.com
Web: <https://www.clinisciences.com>



Denmark

Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: denmark@clinisciences.com
Web: <https://www.clinisciences.com>



Finland

Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: suomi@clinisciences.com
Web: <https://www.clinisciences.com>



France

Company: CliniSciences S.A.S
Address: 74 Rue des Suisses, 92000
Nanterre- France
Telephone: +33 9 77 40 09 09
Fax: +33 9 77 40 10 11
Email: info@clinisciences.com
Web: <https://www.clinisciences.com>



Germany

Company: Clinisciences GmbH
Address: Wilhelm-Mausier-Str. 41-43,
50827 Köln - Germany
Telephone: +49 221 9498 320
Fax: +49 221 9498 325
Email: deutschland@clinisciences.com
Web: <https://www.clinisciences.com>



Iceland

Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: island@clinisciences.com
Web: <https://www.clinisciences.com>



Ireland

Company: CliniSciences Limited
Address: Ground Floor, 71 lower Baggot street
Dublin D02 P593 - Ireland
Telephone: +353 1 6971 146
Fax: +353 1 6971 147
Email: ireland@clinisciences.com
Web: <https://www.clinisciences.com>



Italy

Company: CliniSciences S.r.l
Address: Via Maremmana inferiore 378
Roma 00012 Guidonia Montecelio - Italy
Telephone: +39 06 94 80 56 71
Fax: +39 06 94 80 00 21
Email: italia@clinisciences.com
Web: <https://www.clinisciences.com>



Netherlands

Company: CliniSciences B.V.
Address: Kraaijenhoffstraat 137A,
1018RG Amsterdam, - Netherlands
Telephone: +31 85 2082 351
Fax: +31 85 2082 353
Email: nederland@clinisciences.com
Web: <https://www.clinisciences.com>



Norway

Company: CliniSciences AS
Address: c/o MerVerdi Munkerudtunet 10
1164 Oslo - Norway
Telephone: +47 21 988 882
Email: norge@clinisciences.com
Web: <https://www.clinisciences.com>



Poland

Company: CliniSciences sp.Z.o.o.
Address: ul. Rotmistrza Witolda Pileckiego 67
lok. 200 - 02-781 Warszawa -Poland
Telephone: +48 22 307 0535
Fax: +48 22 307 0532
Email: polska@clinisciences.com
Web: <https://www.clinisciences.com>



Portugal

Company: Quimigen Unipessoal LDA
Address: Rua Almada Negreiros, Lote 5, Loja 14,
2615-275 Alverca Do Ribatejo - Portugal
Telephone: +351 30 8808 050
Fax: +351 30 8808 052
Email: info@quimigen.com
Web: <https://www.quimigen.pt>



Spain

Company: CliniSciences Lab Solutions
Address: C/ Hermanos del Moral 13
(Bajo E), 28019, Madrid - Spain
Telephone: +34 916 750 700
Fax: +34 91 269 40 74
Email: espana@clinisciences.com
Web: <https://www.clinisciences.com>



Sweden

Company: CliniSciences ApS
Address: Oesterbrogade 226, st. 1,
Copenhagen, 2100 - Denmark
Telephone: +45 89 888 349
Fax: +45 89 884 064
Email: sverige@clinisciences.com
Web: <https://www.clinisciences.com>



Switzerland

Company: CliniSciences AG
Address: Fracht Ost Flughafen Kloten
CH-8058 Zürich - Switzerland
Telephone: +41 (044) 805 76 81
Fax: +41 (044) 805 76 75
Email: switzerland@clinisciences.com
Web: <https://www.clinisciences.com>



UK

Company: CliniSciences Limited
Address: 11 Progress Business center, Whittle
Parkway, SL1 6DQ Slough- United Kingdom
Telephone: +44 (0)1753 866 511
or +44 (0) 330 684 0982
Fax: +44 (0)1753 208 899
Email: uk@clinisciences.com
Web: <https://www.clinisciences.com>



USA

Company: CliniSciences LLC
Address: c/o Carr Riggs Ingram,
500 Grand Boulevard, Suite 210 Miramar
Beach, FL 32550- USA
Telephone: +1 850 650 7790
Fax: +1 850 650 4383
Email: usa@clinisciences.com
Web: <https://www.clinisciences.com>

